designed for scientists

LT 5.28

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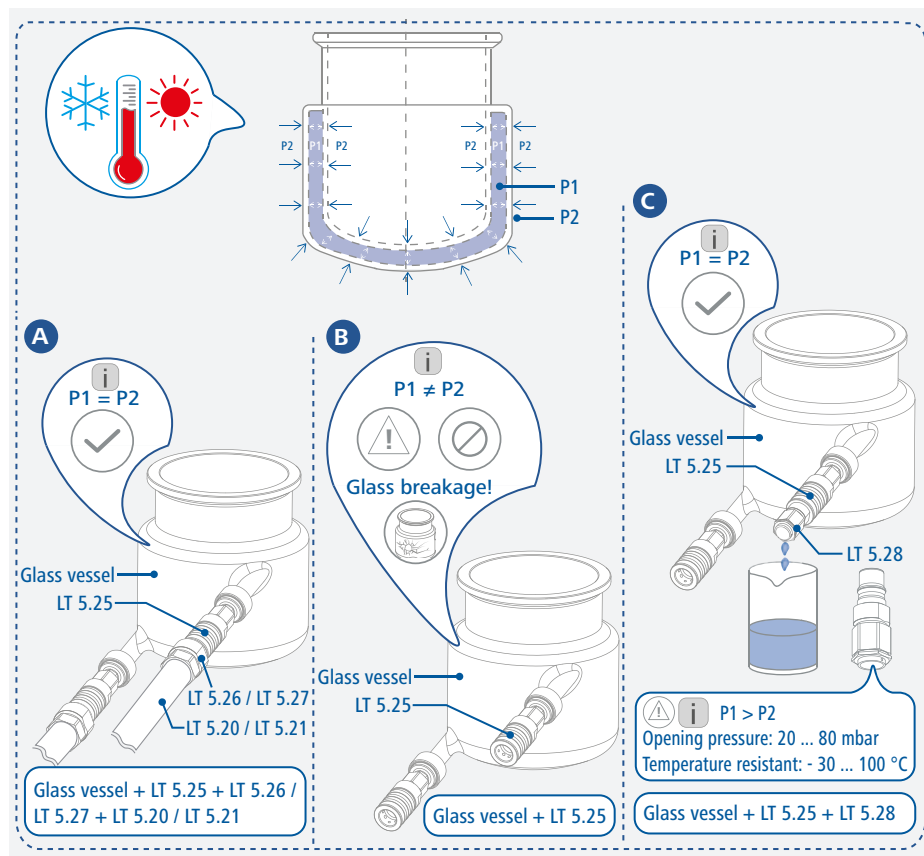
Notice!

- LT 5.28 relief valve serves to prevent glass breakage of the double jacket reaction vessel, especially during assembly and disassembly from the heat transfer fluid circuit and during storage of the filled reaction vessel with LT 5.25 pressure-tight lock coupling assembled.
- The LT 5.28 relief valve is used with the LT 5.25 pressure-tight lock coupling to equalize the pressure in the otherwise hermetically sealed chamber of the reaction vessel with the environment.

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Risk of burns!

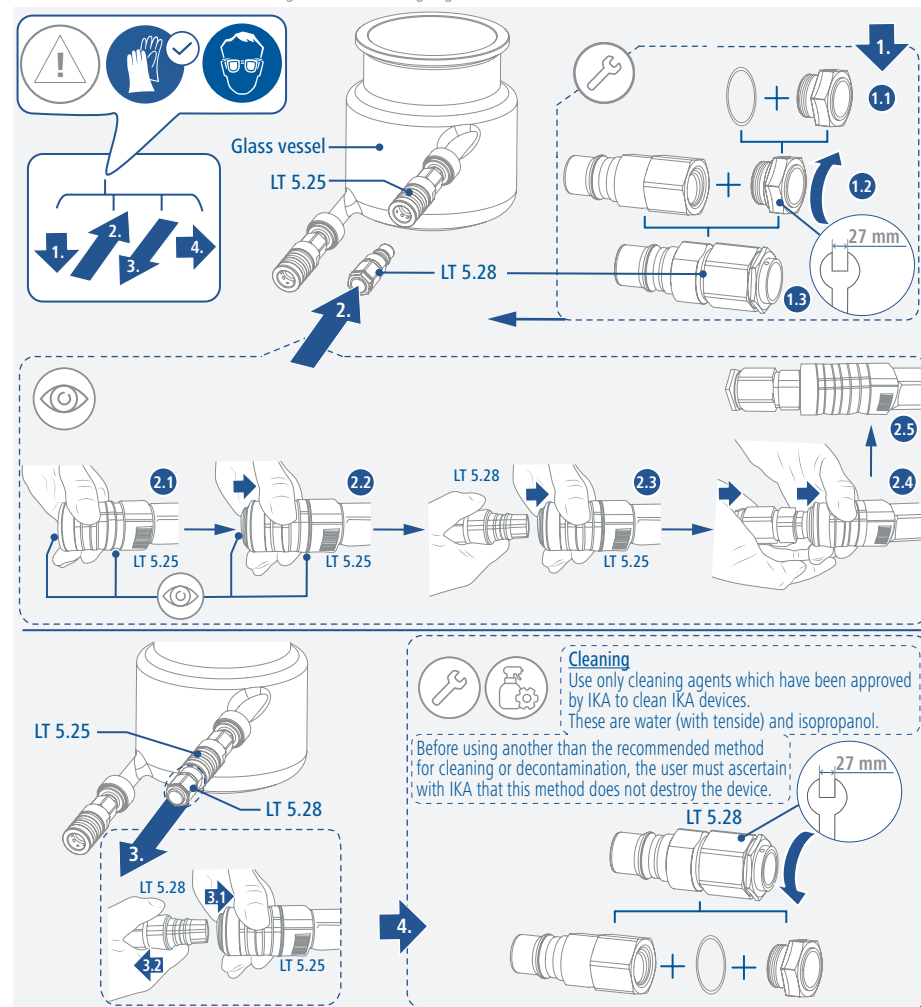
- Reactor must have cooled down to room temperature!

Notice!

- Wear your personal protective equipment during all work!

- Assemble LT 5.28 relief valve.
- Connect LT 5.28 to LT 5.25 lock coupling on the glass vessels. Push and hold LT 5.25 with one hand while pushing LT 5.28 into it with the other hand.
- Remove LT 5.28 from LT 5.25. Push and hold LT 5.25 with one hand while pulling out LT 5.28 with the other hand.
- Disassemble LT 5.28 for thorough cleaning.

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Phone: +84 28 38202142

eMail: sales.lab-vietnam@ika.com

Discover and order the fascinating products of IKA online:
www.ika.com

IKAworldwide

IKAworldwide /// #lookattheblue

@IKAworldwide

Technical specifications may be changed without prior notice.

20000030561_20102591_LT 5.28_012021

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited