IKA-Werke GmbH & Co. KG Janke & Kunkel-Straße 10, 79219 Staufen, Germany Phone: +49 7633 831-0 eMail: sales@ika.de		
USA IKA Works, Inc. Phone: +1 910 452-7059 eMail: sales@ika.net	KOREA IKA Korea Ltd. Phone: +82 2 2136 6800 eMail: sales-lab@ika.kr	BRAZIL IKA Brasil Phone: +55 19 3772 9600 eMail: sales@ika.net.br
MALAYSIA IKA Works (Asia) Sdn Bhd Phone: +60 3 6099-5666 eMail: sales.lab@ika.my	CHINA IKA Works Guangzhou Phone: +86 20 8222 6771 eMail: info@ika.cn	POLAND IKA Poland Sp. z o.o. Phone: +48 22 201 99 79 eMail: sales.poland@ika.com
JAPAN IKA Japan K.K. Phone: +81 6 6730 6781 eMail: info_japan@ika.ne.jp	INDIA IKA India Private Limited Phone: +91 80 26253 900 eMail: info@ika.in	UNITED KINGDOM IKA England LTD. Phone: +44 1865 986 162 eMail: sales.england@ika.com
VIETNAM IKA Vietnam Company Limited Phone: +84 28 38202142 eMail: sales.lab-vietnam@ika.com	THAILAND IKA Works (Thailand) Co. Ltd. Phone: +66 2059 4690 eMail: sales.lab-thailand@ika.com	TURKEY IKA Turkey A.Ş. Phone: +90 216 394 43 43 eMail: sales.turkey@ika.com
Discover and order the fascinating products of IKA online: www.ika.com		
 IKAworldwide	 IKAworldwide /// #lookattheblue	 @IKAworldwide

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/// Warnsymbole

Gefahr!

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Warning symbols

- 

Danger! Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.
- 

Warning! Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.
- 

Caution! Indicates a potentially hazardous situation, which, if not avoided, can result in injury.
- 

Notice! Indicates practices which, if not avoided, can result in equipment damage.

General Symbols

- 

Position number
Indicates device components relevant to actions.
- 

Wrong
Indicates the incorrect execution of an action step.
- 

Correct / result
Indicates the correct execution or the result of an action step.
- 

Note
Indicates steps of actions that require particular attention.

Safety instructions

General information

- ›

Read the operating instructions in its entirety before using the device and follow the safety instructions. If this accessory is used with another device, observe also its operating instructions.
- ›

Keep the operating instructions in a place where it can be accessed by everyone.
- ›

Ensure that only trained staff work with the device.
- ›

Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

Device design

- 

Notice!
- ›

Keep the adhesive surface free of dust.
- ›

Check the adhesive force of the adhesive pad after each removal and replacement.
- ›

Clean the adhesive pad of the stirring attachment with a surfactant-containing water moistened, lint-free cloth.
- ›

Protect the device and accessories from bumps and impacts.
- ›

Check the device and accessories for damage before each use. Do not use damaged components.

Working with the device

- 

Danger!
- ›

To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.
- 

Warning!
- ›

Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- ›

Beware of hazards due to:

-

glass breakage as a result of mechanical shaking power,

-

incorrect container size,

-

overflowing of media,

-

unsafe condition of container.
- ›

When handling hazardous substances or mixtures of substances which are toxic or are contaminated with pathogenic micro-organisms, the user should take appropriate precautions. Observe the national regulations, the biosafety level of your laboratory, the material safety data sheets.
- ›

For germs or biological material or risk group II or higher, please observe the “Laboratory Biosafety Manual” issued by the WorldHealth Organization).

- 

Caution!
- ›

Wear your personal protective equipment in accordance with the hazard category of the media to be processed. There may be a risk from:

-

splashing and evaporation of liquids,

-

ejection of parts,

-

release of toxic or combustible gases.

Accessories

- ›

Safe operation is guaranteed only with the use of original IKA accessories.
- ›

Observe the operating instructions of the accessories.

Maintenance

- ›

The device may only be opened by qualified and IKA approved experts.

Disposal instructions

- ›

The device, accessories and packaging must be disposed of in accordance with local and national regulations.

Intended use

Use

- ›

The TW.IX is a nine-position stirring attachment for mixing of small vessels simultaneously.
- ›

The vessel carriers TW.IX.23 & TW.IX.28 are used for higher safety of stability and convenient and fast exchange of all vessels simultaneously.
- ›

Equip the anti-slip pad TW.IX.ANTISLIP to reduce noise and achieve better self-holding properties of the vessels when stirring with liquids of higher viscosity.

Area of use

- ›

Indoor environments similar to that a laboratory of research, teaching, trade or industry area.
- ›

The safety of the user cannot be guaranteed:

-

if the device is operated with accessories that are not supplied or recommended by the manufacturer,

-

if the device is operated improperly or contrary to the manufacture's specifications,

-

if the device or the printed circuit board are modified by third parties.

Unpacking

Unpacking

- ›

Unpack the device carefully. Any damage should immediately be reported to the carrier (mail, rail or freight forwarding company).

Scope of delivery

TW.IX	
1	TW.IX
2	IKAFLON 15 round (5 pcs.), 2x
4	Operating instructions

TW.IX.23/28	
1	TW.IX.23/28
3	TW.IX.ANTISLIP
4	Operating instructions

Assembly

1.

Degrease the installation surface of the drive unit (e.g. with isopropanol).
2.

Remove the protective film from the adhesive mat.
3.

Place the stirring attachment in the center of the drive unit.
4.

Remove the stirring attachment by rotating it around its own axis. This creates shear forces that detach the attachment from the drive unit.
5.

Place the anti-slip mat at an even distance from the outer edge of the vessel carrier.

Cleaning

- ›

Use only cleaning agents which have been approved by IKA to clean IKA devices. These are: water (containing surfactant) and isopropyl alcohol.
- ›

Wear protective gloves during cleaning the devices.
- ›

Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.
- ›

The adhesive force of the adhesive pad can be restored by cleaning it with water.

Accessories

- ›

For accessories see www.ika.com.

Technical data

TW.IX	
Number of stirring positions	9 pcs
Stirring position distance	30 mm
Stirring quantity min. per stirring position (H ₂ O)	8 ml
Stirring quantity max. per stirring position (H ₂ O)	60 ml
Vessel diameter min.	12 mm
Vessel diameter max.	27.5 mm
Dimensions (W x D x H)	95 x 95 x 21 mm
Weight	0.37 kg

TW.IX.23/28	
Number of stirring positions	9 pcs
Stirring position distance	30 mm
Vessel diameter min.	12 mm
Vessel diameter max.	23/27.5 mm
Dimensions (W x D x H)	99 x 101 x 47.5 mm
Weight	0.2 kg

- ›

Subject to technical changes!