



designed for scientists


IKA ICC basic IB R RO 15 eco
IKA ICC control IB R RO 15 eco




Observe the Operating Instructions of the ICC basic / control and RO 15!




Delivery scope

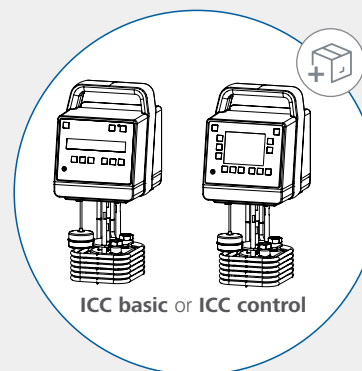


ICC basic or ICC control

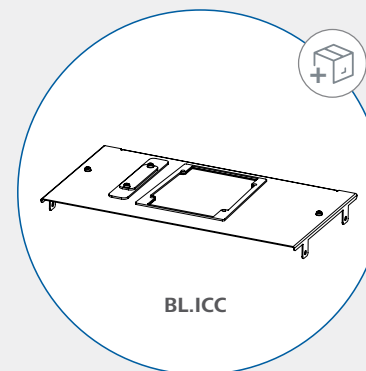


BL.ICC

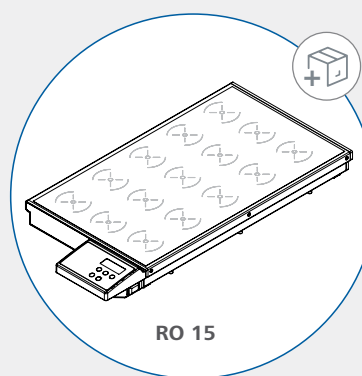


RO 15

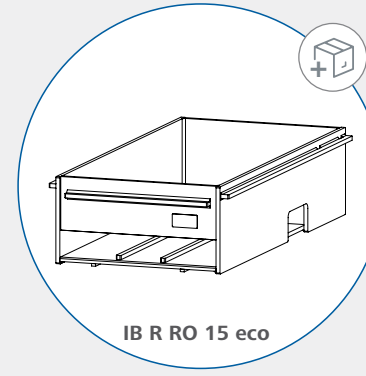

IB R RO 15 eco




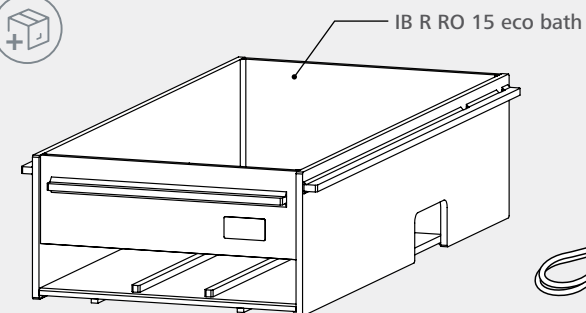



Extension cord

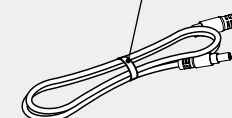







Installation

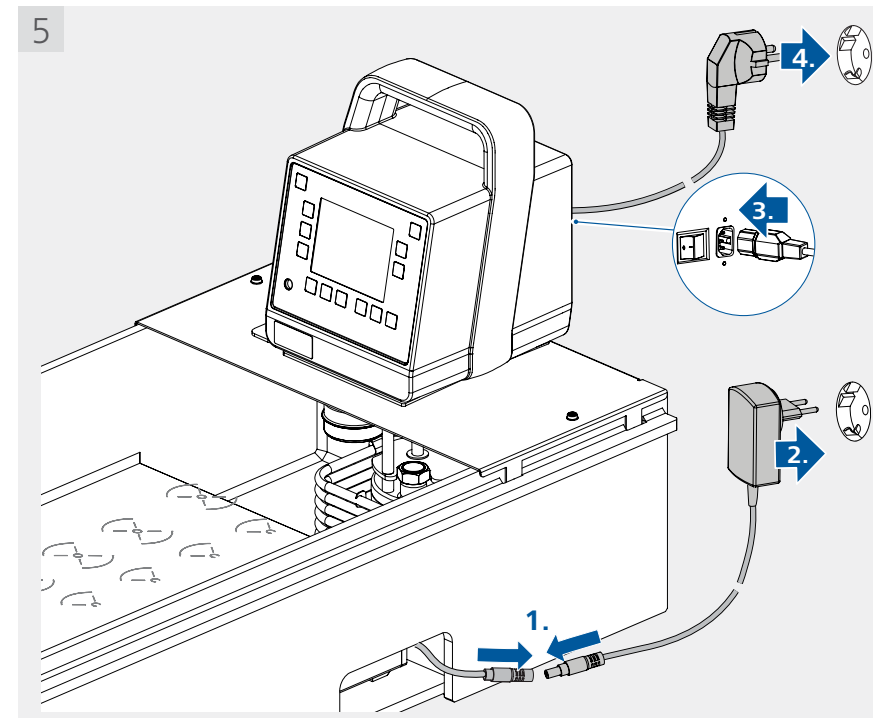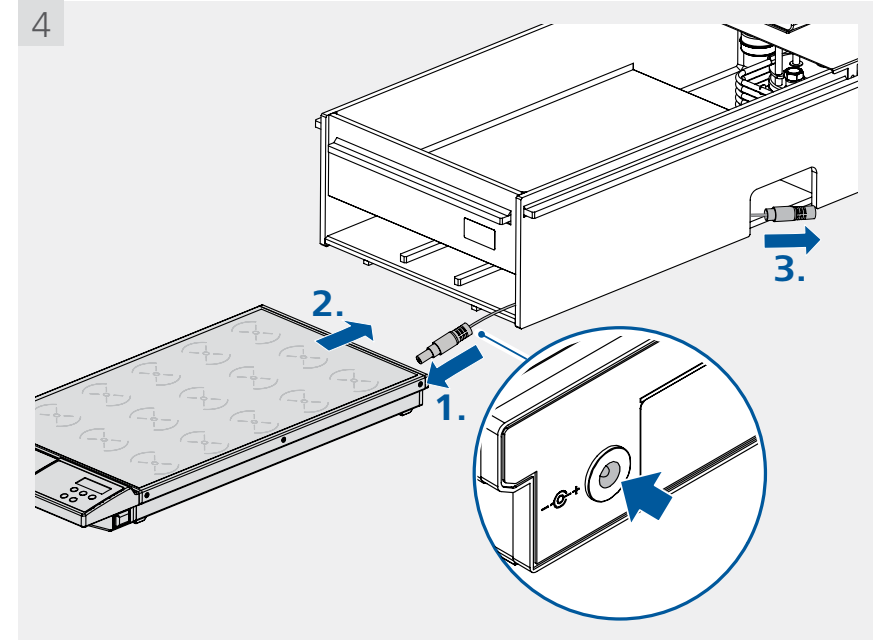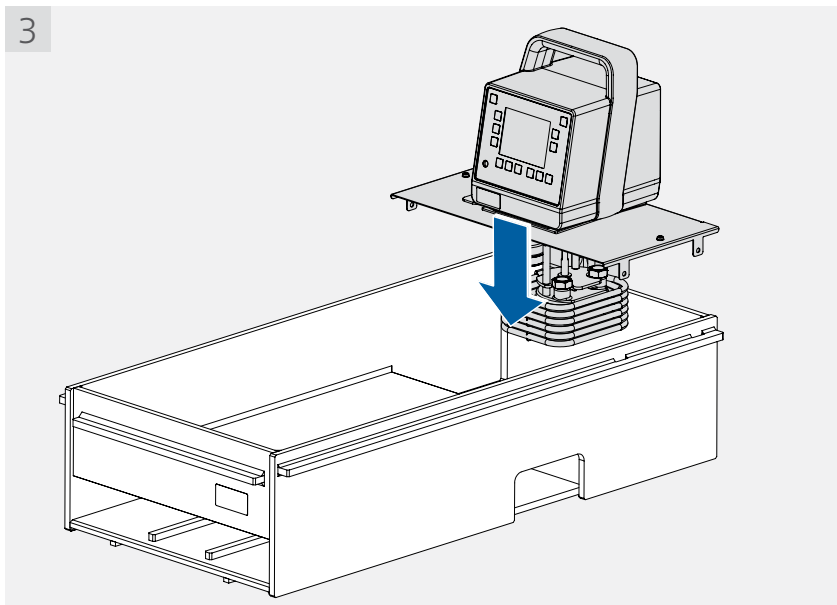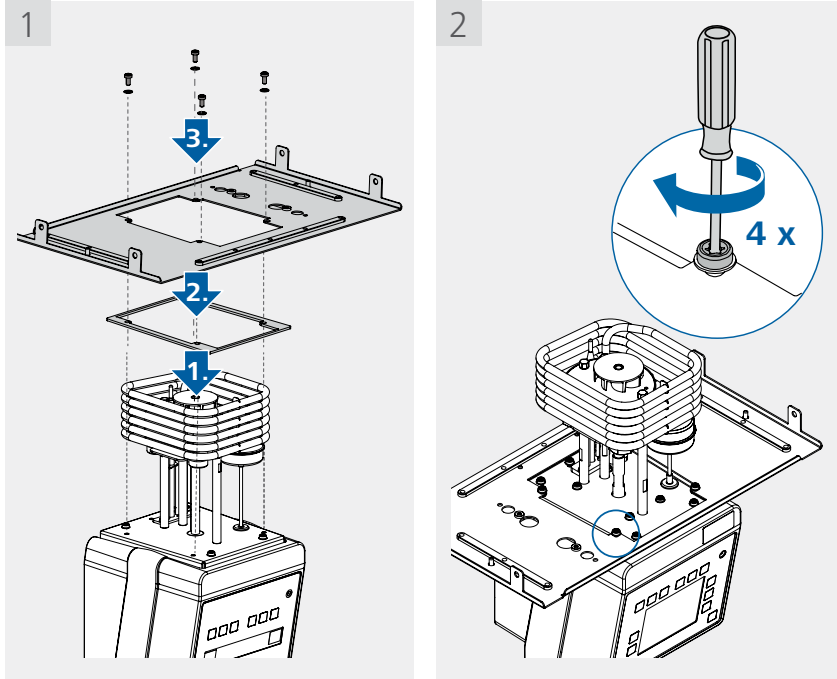

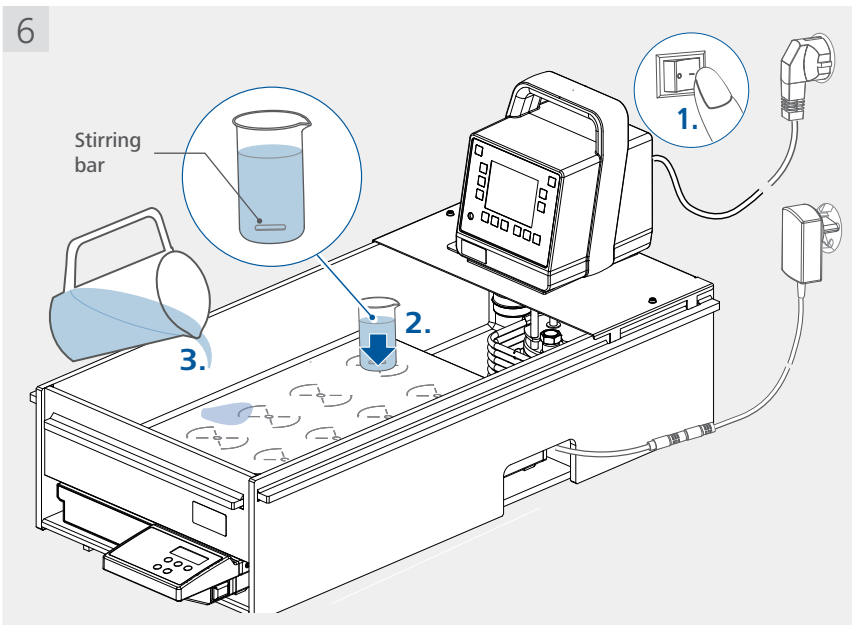


! Untreated tap water is not recommended. It is recommended to use distilled water or high purity water (ion exchangers) and add 0.1 g soda (sodium carbonate Na_2CO_3) / liter, to reduce corrosive properties.

i In order to improve the temperature uniformity of the fluid in the bath, you can connect the accessory pump distribution set **PDS.IB.R.RO**.

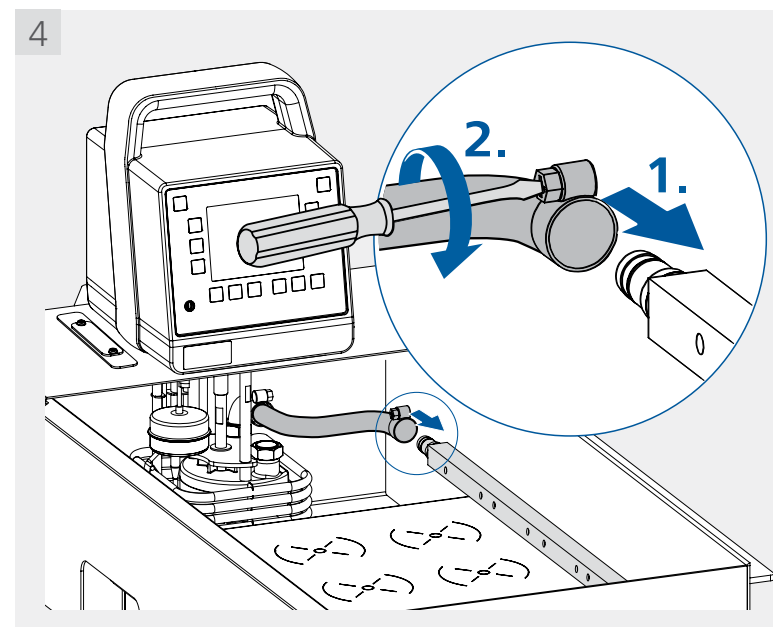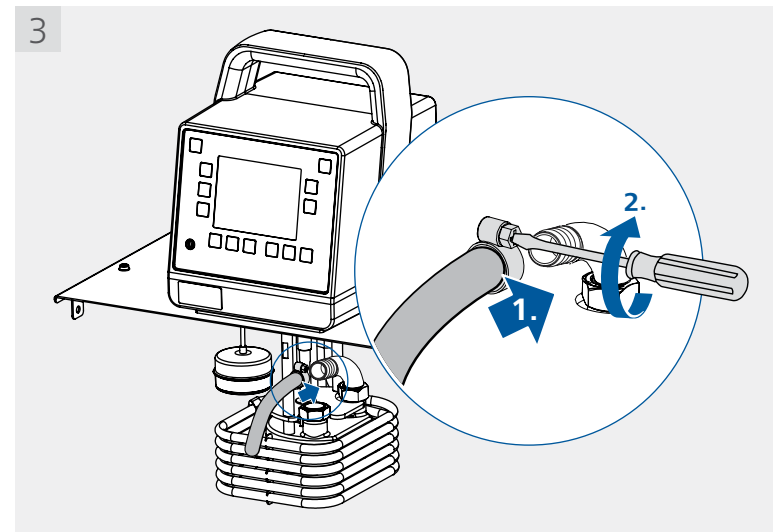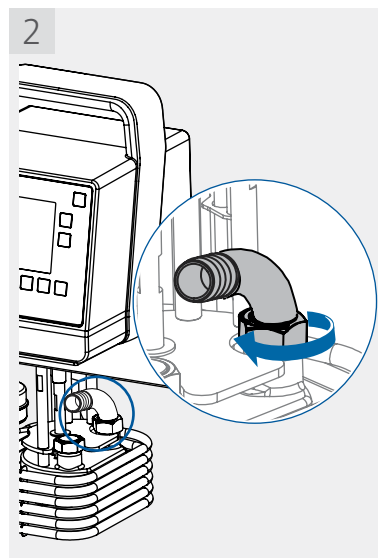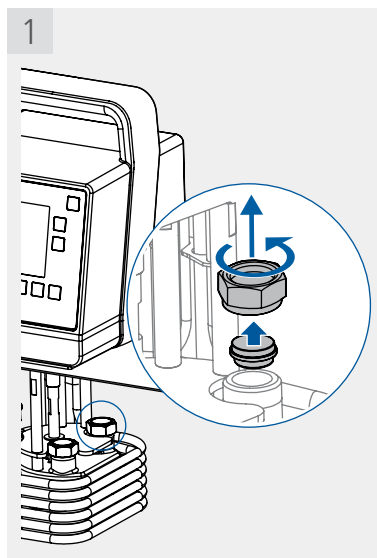



Operation

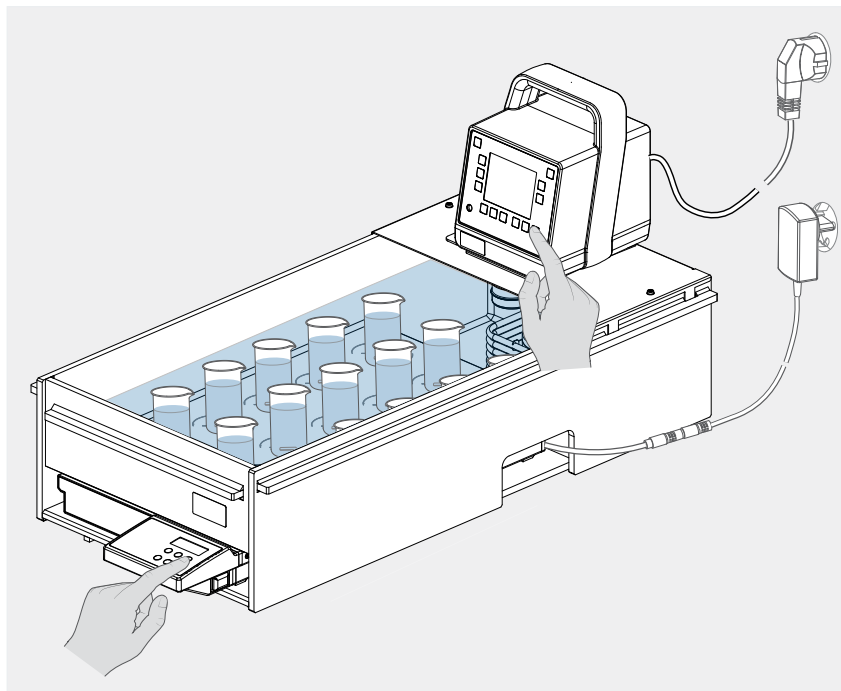

For operation, please refer to the operating instructions of the **ICC basic / control** and **RO 15**!




The maximum temperature in the plastic bath (**IB R RO15 eco**): 80 °C (water only)!


For higher temperature and oil use, we recommend the stainless steel bath (**IB R RO15 pro**) (max. 130 °C).

Note: when working at temperatures higher than 45 °C, we suggest to use floating globes to avoid too fast loss of water by evaporation. Maximum temperature difference between refilled water and actual bath temperature should be less than 10 K.

Transportation:

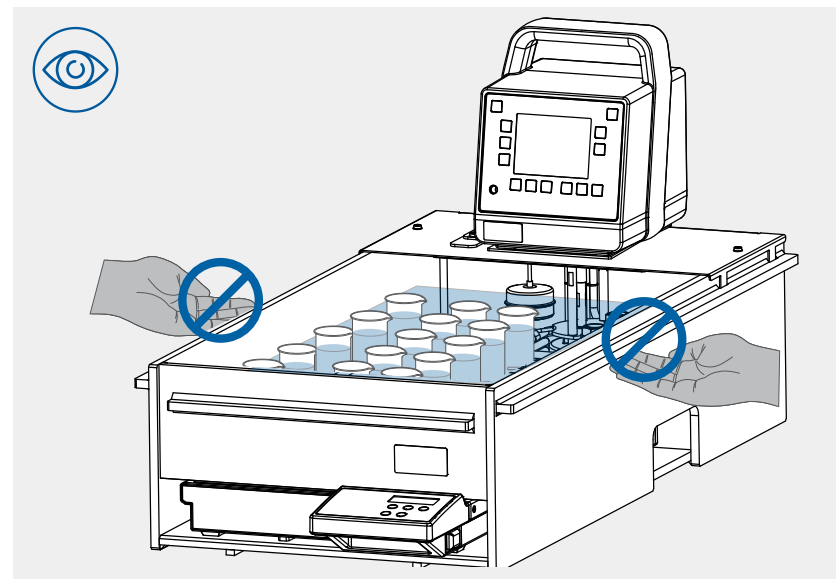

Before transportation, please remove the devices (**ICC & RO 5**) and empty the fluid in the bath! Never lift the handles when the bath is filled with fluid!

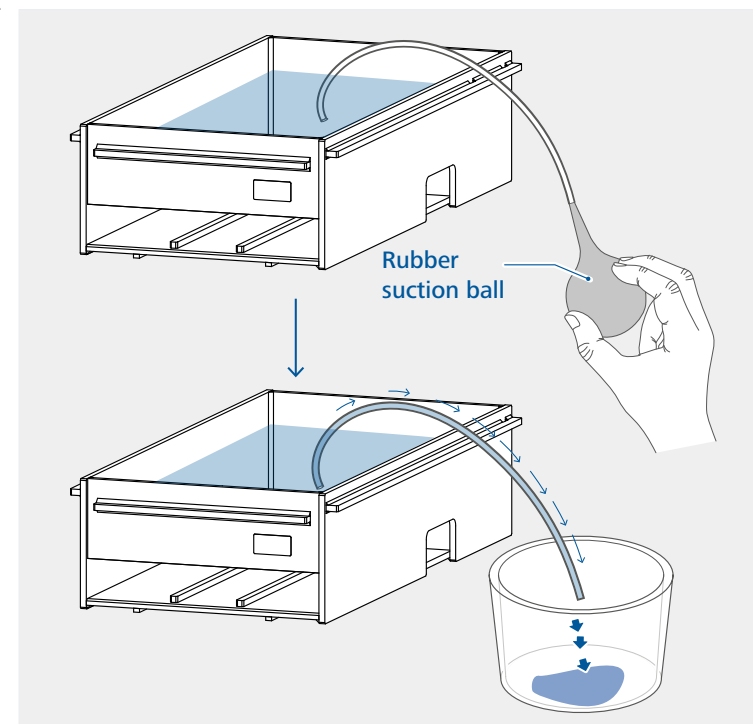

Note:



designed for scientists

IKA-Werke GmbH & Co. KG

Janke & Kunkel-Straße 10, 79219 Staufen, Germany
Phone: +49 7633 831-0, Fax: +49 7633 831-98
eMail: sales@ika.de

USA

IKA Works, Inc.
Phone: +1 910 452-7059
sales@ika.net

KOREA

IKA Korea Ltd.
Phone: +82 2 2136 6800
sales-lab@ika.kr

BRASIL

IKA Brasil
Phone: +55 19 3772-9600
sales@ika.net.br

MALAYSIA

IKA Works (Asia) Sdn Bhd
Phone: +60 3 6099-5666
sales.lab@ika.my

CHINA

IKA Works Guangzhou
Phone: +86 20 8222 6771
info@ika.cn

POLAND

IKA Poland Sp. z o.o.
Phone: +48 22 201 99 79
sales.poland@ika.com

JAPAN

IKA Japan K.K.
Phone: +81 6 6730 6781
info_japan@ika.ne.jp

INDIA

IKA India Private Limited
Phone: +91 80 26253 900
info@ika.in

UNITED KINGDOM

IKA England LTD.
Phone: +44 1865 986 162
sales.england@ika.com

VIETNAM

IKA Vietnam Company Limited
Phone: +84 28 38202142
sales.lab-vietnam@ika.com

THAILAND

IKA Works (Thailand) Co. Ltd.
Phone: +66 2059 4690
sales.lab-thailand@ika.com

TURKEY

IKA Turkey A.S
Phone: +90 216 394 43 43
sales.turkey@ika.com

KENYA

IKA Works Kenya Ltd.
Phone: +254 112 323 745
sales.kenya@ika.com

UGANDA

IKA Works Kampala Limited
Phone: +254 112 323 745
sales.uganda@ika.com

SPAIN

IKA Works Spain, S. L.
Barcelona
sales.spain@ika.com

Discover and order the fascinating products of IKA online:
www.ika.com



IKAWorldwide



IKAWorldwide /// #lookattheblue



@IKAWorldwide

Technical specifications may be changed without prior notice.