IKA-Werke GmbH & Co. KG
Janke & Kunkel-Straße 10, 79219 Staufen, Germany
Phone: +49 7633 831-0
eMail: sales@ika.de

USA
IKA Works, Inc.
Phone: +1 910 452-7059
eMail: sales@ika.net

MALAYSIA
IKA Works (Asia) Sdn Bhd
Phone: +60 3 6099-5666
eMail: sales.lab@ika.my

JAPAN
IKA Japan K.K.
Phone: +81 6 6730 6781
eMail: info_japan@ika.ne.jp

VIETNAM
IKA Vietnam Company Limited
Phone: +84 28 38202142
eMail: sales.lab-vietnam@ika.com

KENYA
IKA Works Kenya Ltd.
Phone: +254 112 323 745
eMail: sales.kenya@ika.com

KOREA
IKA Korea Ltd.
Phone: +82 2 2136 6800
eMail: sales-lab@ika.kr

CHINA
IKA Works Guangzhou
Phone: +86 20 8222 6771
eMail: info@ika.cn

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Explication of warning symbols

Warning symbols

- **Danger!** Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.
- **Warning!** Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.
- **Caution!** Indicates a potentially hazardous situation, which, if not avoided, can result in injury.
- **Notice!** Indicates practices which, if not avoided, can result in equipment damage.

General Symbols

- **A** Position number
Indicates device components relevant to actions.
- **Wrong**
Indicates the incorrect execution of an action step.
- **Correct / result**
Indicates the correct execution or the result of an action step.
- **Note**
Indicates steps of actions that require particular attention.

Safety instructions

General information

- Read the operating instructions in its entirety before using the device and follow the safety instructions. If this accessory is used with another device, observe also its operating instructions.
- Keep the operating instructions in a place where it can be accessed by everyone.
- Ensure that only trained staff work with the device.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

Device design

- **Notice!**
- Keep the adhesive surface free of dust.
 - Check the adhesive force of the adhesive pad after each removal and replacement.
 - Clean the adhesive pad of the shaking attachment with a surfactant-containing water moistened, lint-free cloth.
 - Protect the device and accessories from bumps and impacts.
 - Check the device and accessories for damage before each use. Do not use damaged components.

Working with the device

- **Danger!**
- To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.
- **Warning!**
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
 - Beware of hazards due to:
 - flammable materials,
 - glass breakage as a result of mechanical shaking power, incorrect container size,
 - overflowing of media,
 - unsafe condition of container.
 - Keep hand/arm vibration to a minimum (Directive 2002/44/EC) when operated manually:
 - Keep the sample container as vertical as possible.
 - Only exert the minimum pressure necessary for the mixing process.
 - Use attachments for multiple samples when there is a large number of samples.
 - If the shaking movement of the appliance increases (resonance), reduce the speed or pass through the critical phase as quickly as possible.
 - Reduce the speed if:
 - the medium splashes out of the vessel because the speed is too high,
 - the appliance is not running smoothly,
 - the appliance begins to move around because of dynamic forces.

- Firmly secure the accessories and vessels in place, otherwise shaking vessels could be damaged or projected out.
- Position one single shaking vessel in the centre and several shaking vessels so that they are evenly spread out.
- Sharp edged vessels cause wear debris on the attachments.
- When handling hazardous substances or mixtures of substances which are toxic or are contaminated with pathogenic micro-organisms, the user should take appropriate precautions. Observe the national regulations, the biosafety level of your laboratory, the material safety data sheets.
- For germs or biological material or risk group II or higher, please observe the “Laboratory Biosafety Manual” issued by the WorldHealth Organization).

- **Caution!**
- Wear your personal protective equipment in accordance with the hazard category of the media to be processed. There may be a risk from:
 - splashing and evaporation of liquids,
 - ejection of parts,
 - release of toxic or combustible gases.

Accessories

- Safe operation is guaranteed only with the use of original IKA accessories.
- Observe the operating instructions of the accessories.

Disposal instructions

- The device, accessories and packaging must be disposed of in accordance with local and national regulations.

Intended use

Use

- The TW.VX is a shaking attachment for mixing liquids.
 - Touch mode for single test tubes.
- Continuous mode for single or several vessels.

Area of use

- Indoor environments similar to that a laboratory of research, teaching, trade or industry area.
- The safety of the user cannot be guaranteed:
 - if the device is operated with accessories that are not supplied or recommended by the manufacturer,
 - if the device is operated improperly or contrary to the manufacture’s specifications,
 - if the device or the printed circuit board are modified by third parties.

Unpacking

Unpacking

- Unpack the device carefully. Any damage should immediately be reported to the carrier (mail, rail or freight forwarding company).

Scope of delivery

1	TW.VX	4	MS 1.21 One-hand insert
2	MS 3.1 Standard attachment	5	Lever tool
3	MS 3.3 Universal attachment	6	Operating instructions

Assembly

- Degrease the installation surface of the drive unit (e.g. with isopropanol).
- Remove the protective film from the adhesive mat.
- Place the shaking attachment in the center of the IKA TWISTER drive unit.
- Remove the shaking attachment with the lever tool, by lifting it upwards.

Cleaning

- Use only cleaning agents which have been approved by IKA to clean IKA devices. These are: water (containing surfactant) and isopropyl alcohol.
- Wear protective gloves during cleaning the devices.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.
- The adhesive force of the adhesive pad can be restored by cleaning it with water.

Accessories

- Permitted speed ranges of the attachments:

Speed range (rpm)		
0/100 ... 2500	0/100 ... 1300	0/100 ... 500
MS 3.1	MS 3.3, MS 3.5, MS 3.51, MS 1.21, MS 1.30, MS 1.31, MS 1.32, MS 1.33, MS 1.34, MS 3.4	STICKMAX II




- For accessories see www.ika.com.

Technical data

Type of movement	orbital
Shaking stroke	4.5 mm
Permissible shaking weight (incl. attachment)	0.5 kg

Maintenance

- The device may only be opened by qualified and IKA approved experts.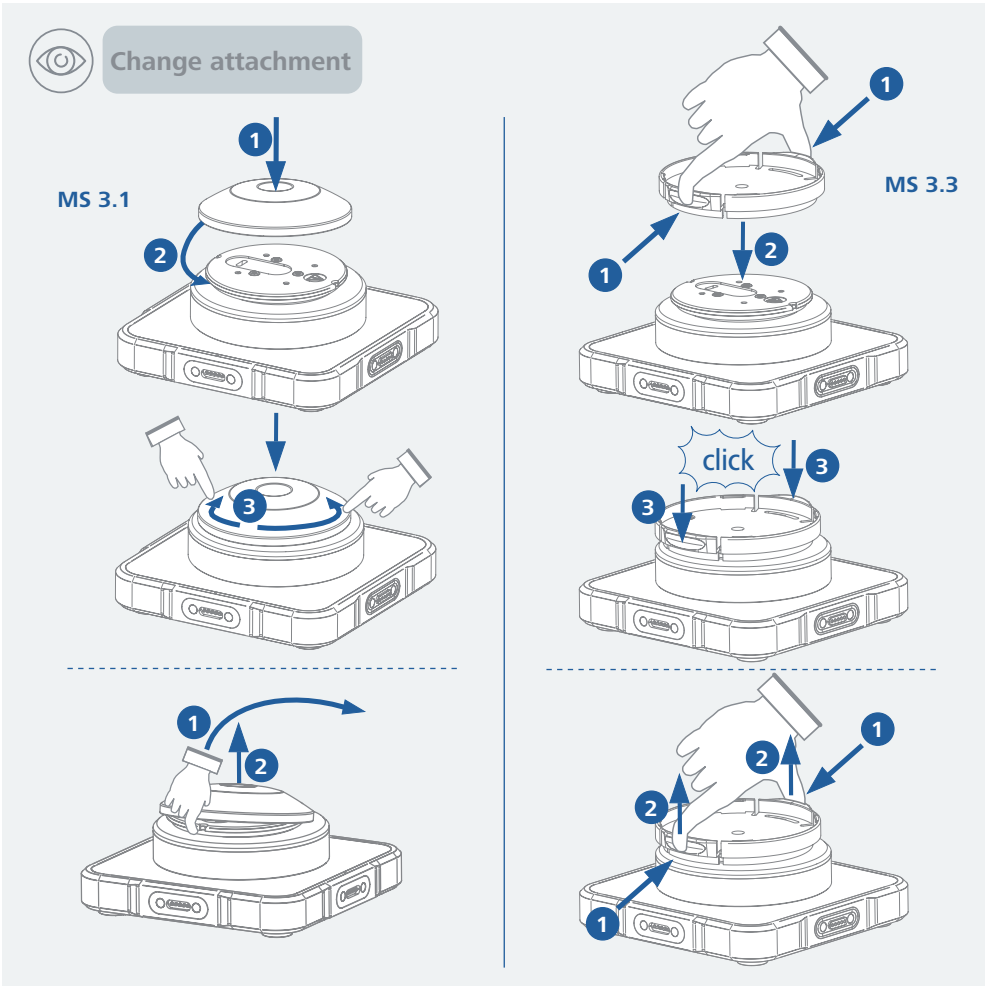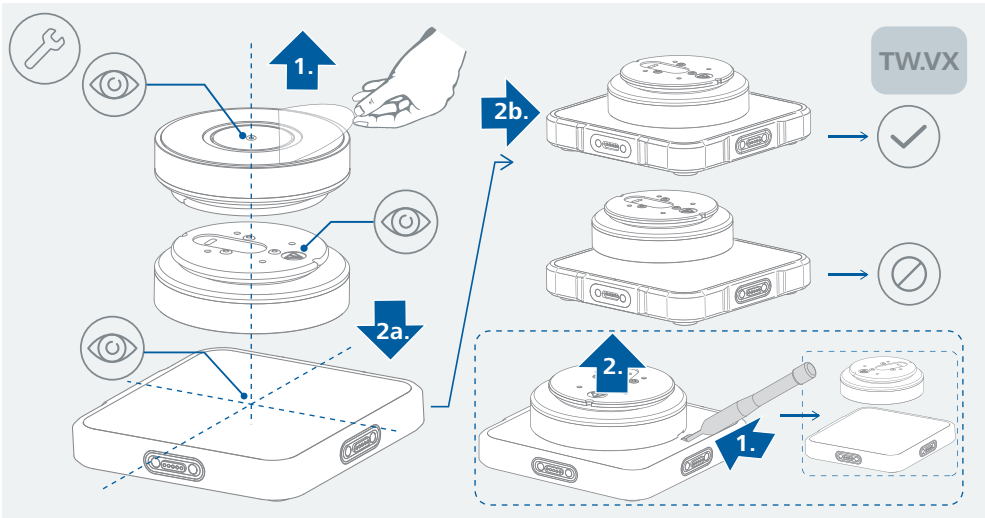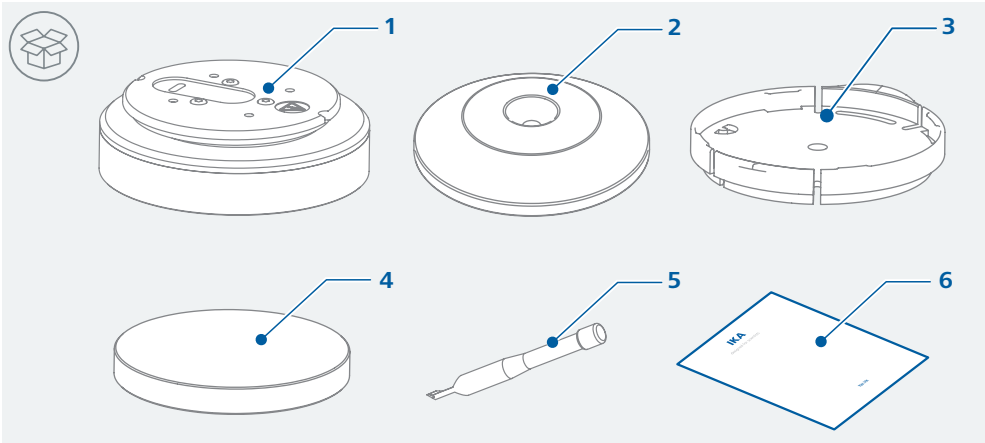

› Subject to technical changes!