designed for scientists

HA.ls.2
HA.ls.10

5
/// DEUTSCH

13
/// FRANÇAIS

21
/// ITALIANO

29
/// POLSKI

37
/// TÜRKÇE

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9
/// ENGLISH

17
/// ESPAÑOL

25
/// PORTUGUÊS

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HA.Is.2

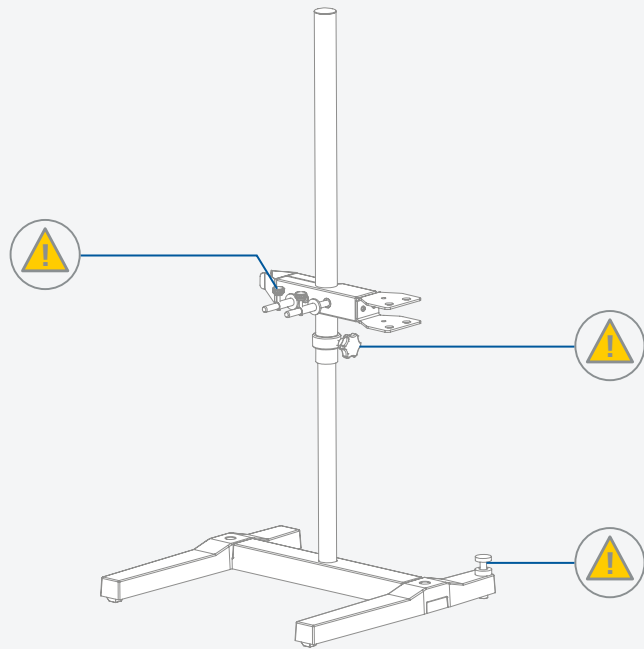
Fig. 1

HA.Is.10

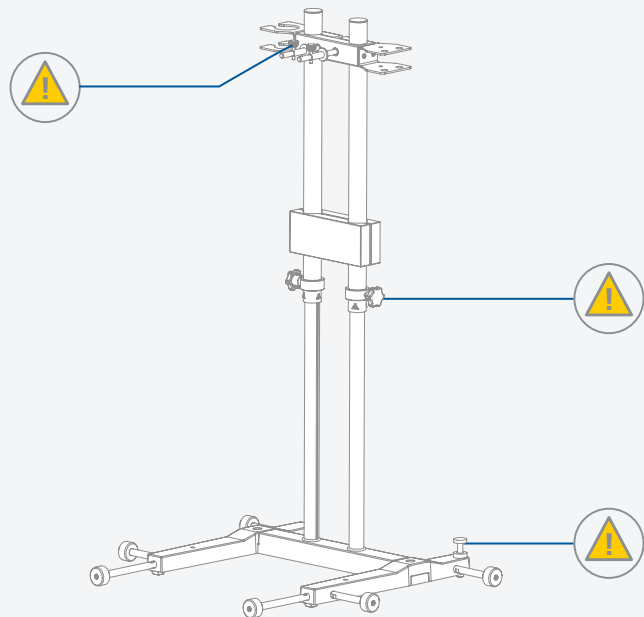
Fig. 2

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	EU Declaration of conformity	9
	Explication of warning symbols	10
	Safety instructions.....	10
	Intended use.....	11
	Unpacking	48
	Assembly	49
	Cleaning	51
	Technical Data	51

Source language: German

EU Declaration of conformity

We declare under our sole responsibility that the product to which this declaration relates is in conformity with directive 2006/42/EC and the following standard or other normative document: EN ISO 12100.

A copy of the complete Declaration of Conformity or further declarations of conformity can be requested at sales@ika.com.

Explication of warning symbols

/// Warning symbols

Danger!

Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.

Warning!

Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.

Caution!

Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Notice!

Indicates practices which, if not avoided, can result in equipment damage.

/// General Symbols

A — Position number

Correct / Result

Shows the correct execution or the result of an action step.

Wrong

Shows the incorrect execution of an action step.

Note

Displays action steps that require particular attention to detail.

Closed lock

Indicates the status of "locked".

Open lock

Indicates the status of "Unlocked".

- › You must also follow the safety instructions contained in the operating instructions for the Bioreactors IKA HABITAT cell/ferment and all other variants.
- › Before use (at regular intervals), check the positioning of the IKA Lid stand to the IKA Habitat Vessel holder (Fig. 7, "Assembly").

Caution!

- › Check the mounting position of the stirrer / disperser to avoid the danger of the unit tipping over.
- › Observe the maximum permissible load-bearing capacity (see "Technical data").
- › The adjustment screw can be used to compensate for unevenness on the set-up plate. Use the lock nut to lock the adjustment screw.
- › When adjusting the height of the movable stand rod, ensure that nothing becomes caught between the devices or other objects attached to the bracket, e.g. the lower end stop or lip of the container.
- › The stand rod must be held in place when loosening the knurled screw (avoid rapid increases).
- › Always store the stand in a relaxed state.

Notice!

- › Pay attention to the marked sites in Fig. 1 and Fig. 2.

Intended use

- › The HA.Is.2 and the HA.Is10 are used only for holding, raising and lowering IKA reactor lids of the IKA HABITAT series.

Notice!

- › The HA.Is.2 may only be used with the IKA reactor lid size 2 l.
- › The HA.Is.10 may only be used with the IKA reactor lid size 5 l and 10 l.
- › The safety of the user cannot be ensured:
 - if the stand is used in conjunctions with accessories not made or recommended by the manufacturer
 - if the stand is not used for the intended purpose as specified by the manufacturer.

Safety instructions

/// General information

- › **Read the operating instructions completely before starting up and follow the safety instructions.**
- › Keep the operating instructions in a place where it can be accessed by everyone.
- › Ensure that only trained staff work with the device.
- › Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- › Install the stand on a flat, stable, clean, non-slip, dry and fire-resistant surface.
- › The device must only be used in a technically perfect condition. The correct functioning of the stand must be checked at regular intervals.
- › Prior to each use, always check the stand for damage. Never use damaged parts.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited
IKA Works (Thailand) Co. Ltd.