
IKA

designed for scientists

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Cleaning tool IKA TS 25-3

Outil de nettoyage IKA TS 25-3

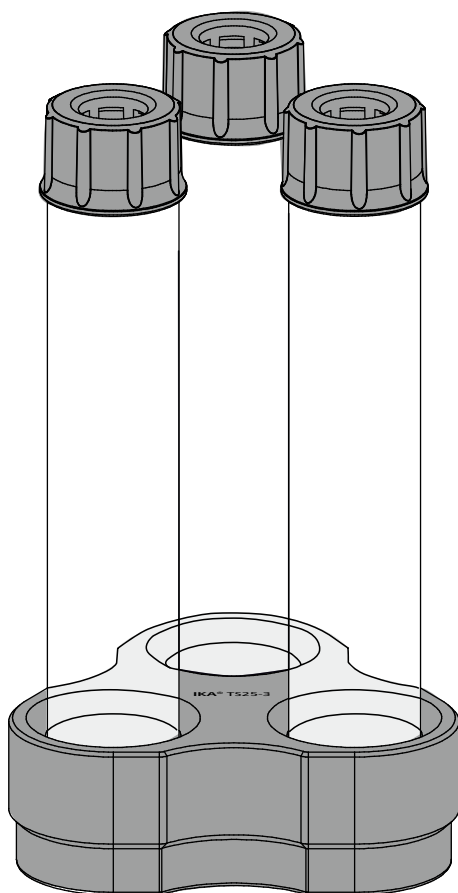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DE

4

Operating instructions





EN

6

Mode d'emploi

FR

8

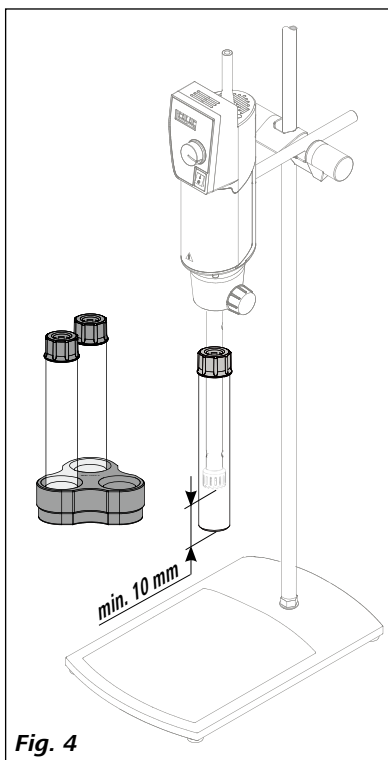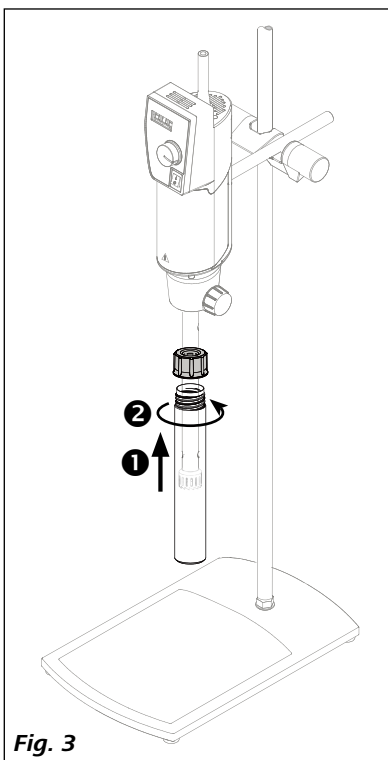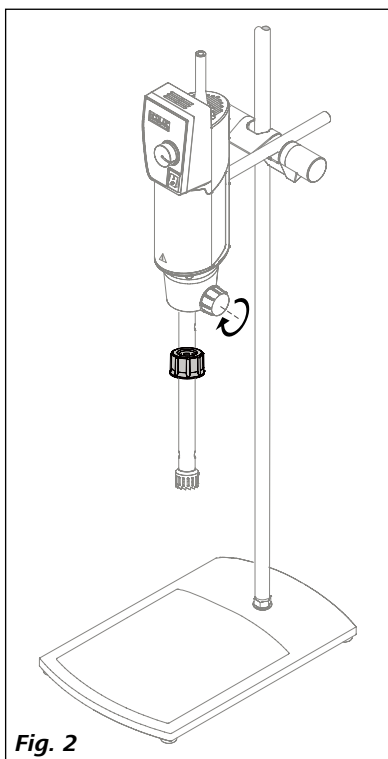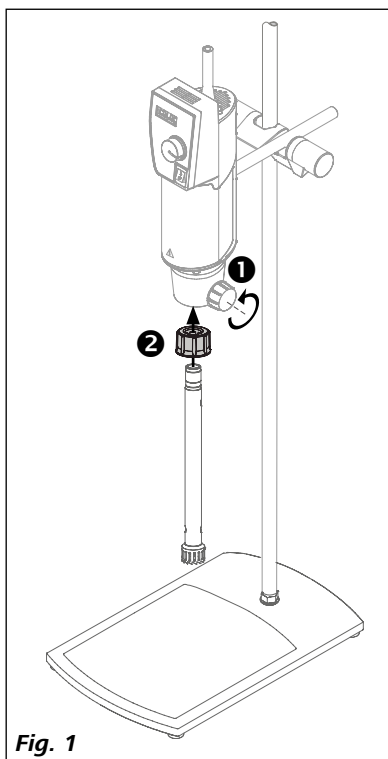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Content

	Page
Explication of warning symbols	6
Safety instructions	6
Correct use	7
Unpacking	7
Product information	7
Installing the cleaning tube to disperser	7
Accessories	7
Technical data	7

Explication of warning symbols

Indicates an (extreme) hazardous situation, which, if not avoided, will result in death, serious injury.

Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.

Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Indicates practices which, if not avoided, can result in equipment damage.

Safety instructions

• Read the operating instructions in full before starting up and follow the safety instructions.

- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the cleaning tube.
- Follow the safety instructions, guidelines, occupational health, safety and accident prevention regulations.
- Set up the tube holder on an even, stable, clean, non-slip, dry and fireproof surface.
- Note the operating instructions of the disperser.

Never run dispersion tools without enough cleaning liquid or medium in the tubes, as the gasket and bearings will be destroyed if the tools are not cooled by the medium.

- Do not process any flammable or combustible materials with the cleaning tubes.
- Only process media that will not react dangerously to the extra energy produced through processing with the tubes. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- A cleaning/dispersing time of a few minutes is usually sufficient. Long processing time will increase the temperature of the cleaning liquid or the medium.
- The tubes are not sealed. The filled tubes are not allowed to lay down!

Wear your personal protective equipment in accordance with the hazard category of the media to be processed. Otherwise there is a risk from:

- splashing and evaporation of liquids
- body parts, hair, clothing and jewellery getting caught.
- Make sure that the stator has always minimum 10mm distance to the tube bottom (see **Fig. 4**).
- The liquid in the tube can get very hot after a short operation. Handle the tube with your personal protective equipment.

Correct use

- **Use:**

Use **TS 25-3** for cleaning of dispersing elements and dispersing of low quantum.

- **Range of use:**

- Laboratories - Schools - Pharmacies - Universities

Unpacking

- **Unpacking:**

- Please unpack the device carefully.
- In the case of any damage a fact report must be sent immediately (post, rail or forwarder).

- **Delivery scope:**

- **TS 25-3** (one tube holder, three tubes and three caps).
- Operating instructions.

Product information

- When use **TS 25-3** in combination with **IKA® T 25 digital** disperser, it can be for cleaning dispersing elements or for dispersing low quantity liquid media in batches.

When use **TS 25-3** for cleaning dispersing elements, we recommend to pre clean the dispersing tool in one tube with suitable cleaning liquid. Use the second tube to wash up the dispersing tool. The third tube is for the final cleaning. The tube holder can be used for holding the tubes.

- **Materials:**

Tubes: PP **Tube cap:** PP **Tube holder:** Aluminum

Installing the cleaning tube to disperser

For installing the tube to disperser, see **Fig. 1** to **Fig. 4**.

Refer to the operating instructions of the disperser.

CAUTION

The tubes are not sealed. Please don't lay them down if cleaning liquid or medium inside.

Accessories

CT.15 15 cleaning tubes with caps

Technical data

Max temperature of the liquid:	70 °C
Min working volume:	50 ml
Max working volume:	100 ml
Suitable for dispersing tools:	all S 25 dispersing tools with Ø 18mm shaft tube

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IKA Korea Ltd.
IKA India Private Limited
IKA England LTD.